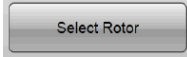
Avanti JXN Series

High-Performance Centrifuge Instructions

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AVANTI JXN SERIES CENTRIFUGES OPERATION

OPERATION STEPS

1. Turn on the power switch on the right side panel.
(If Door wasn't opened, step on the foot pedal to open the door.)
2. Installation Rotor and securely attach the rotor lid knob to the drive shaft by turning it to the right (clockwise). Close the centrifuge door firmly.
3. Start on the Home Page.
 - a. Set the Speed and Rotor :
Select the **SET SPEED** Display/Button. When the Set Speed Page appears, Select the **SELECT ROTOR** Button first to go to the Select Rotor page. When the rotor is selected, Press the **OK** Button to return to the Set Speed Page to set the speed.
 - b. Set the Time :
Select the **SET TIME** Display/Button.
 - i. The maximum time of 99 hours and 59 minutes has been reached.
 - ii. In **HOLD** mode, the run does not end until you select the **STOP** key.
 - iii. The **DELAY START** button to set a future start or stop time. Select **START AT** or **STOP AT** to set a starting or stopping time. Then you can set the date and time in the fields above.
 - c. Set the Temperature :
Select the **SET TEMP** Display/Button, from -10 to $+40^{\circ}\text{C}$.
 - d. Set the Acceleration and Deceleration :
Select the **ACCEL** and **DECEL** Display/Button to select the desired Acceleration and Deceleration profiles. The acceleration value of zero is the maximum (no reduction). The acceleration value of 10 is the slowest (maximum reduction). The same applies to the deceleration values. The value 11 is an absolute reduction, eliminating all braking and allowing the rotor to coast to a stop.
4. Check that all parameters are correct and that the door is closed. Press **START** starting a run and **WAITING TO SET SPEED IS REACHED**.
5. Wait for the set time to count down to zero or end the run by pressing **STOP**. The run is complete, foot pedal to open the door and unmount the rotor and perform all the cleanup.

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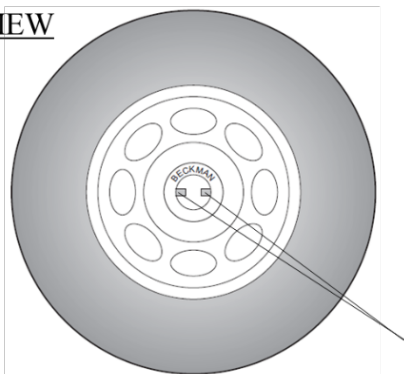
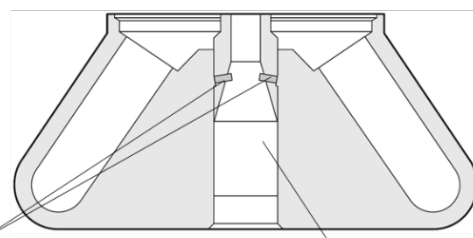
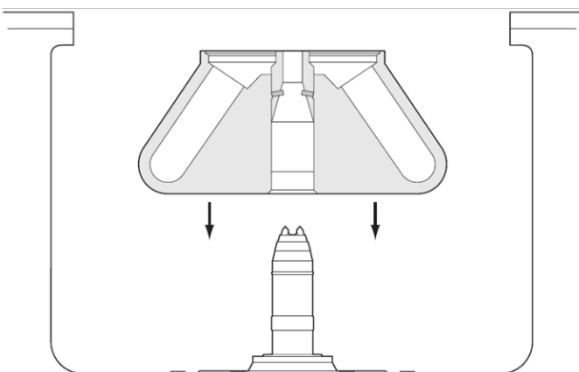
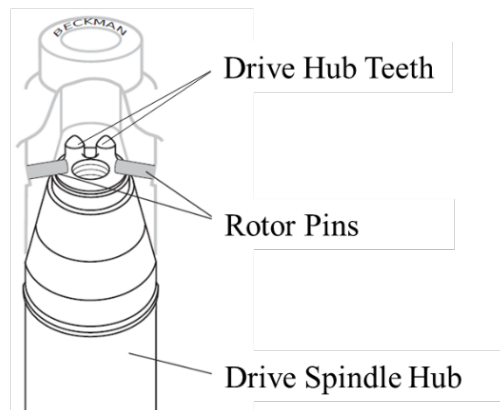
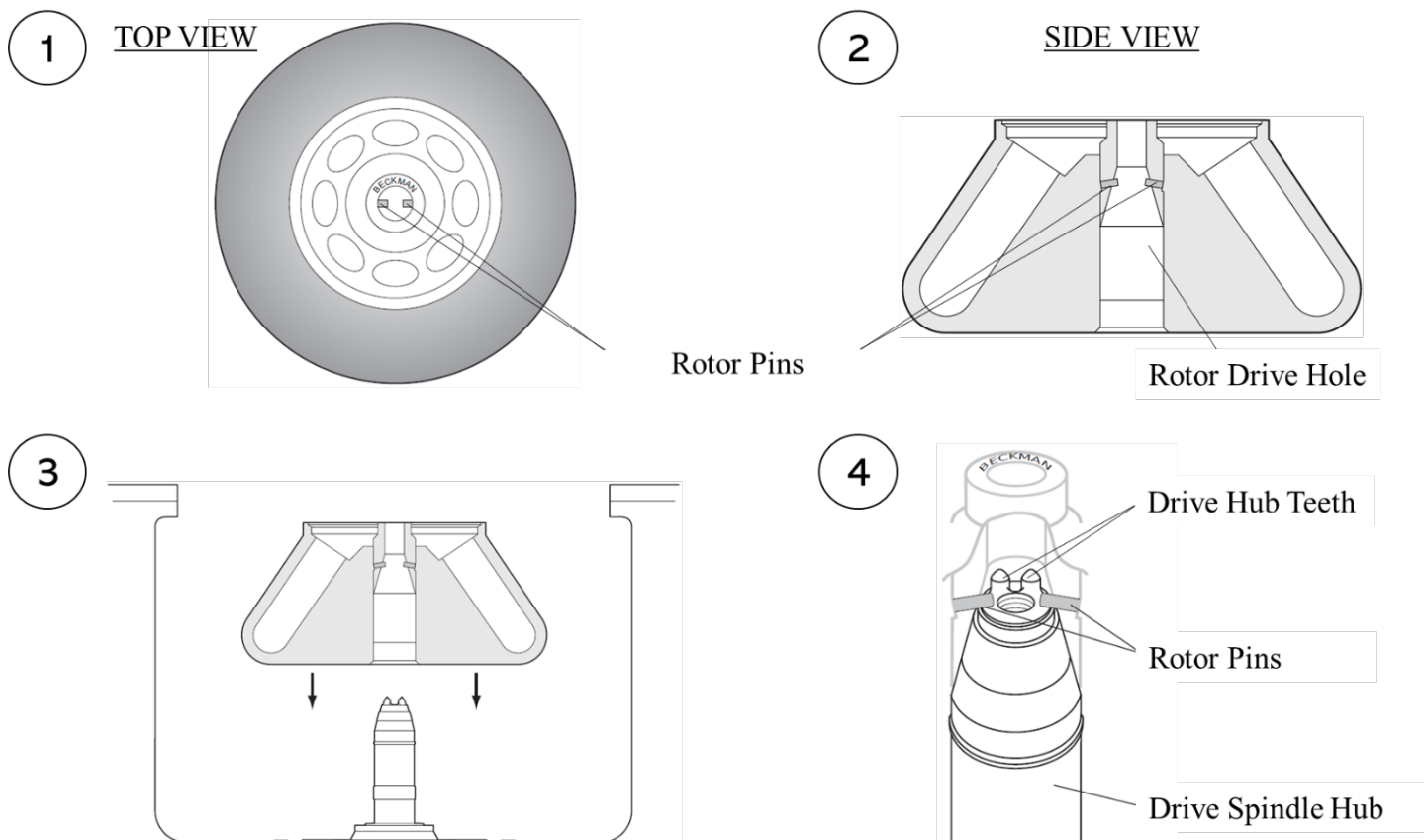
AVANTI J SERIES ROTOR OPERATION

1. Before installing the rotor, please check:
 - a. Inspect the rotor cavities, for rough spots or pitting, white powder deposits. If any of these signs are evident, wash the rotor and rotor components immediately.
 - b. Make sure the rotor and lid's o-rings are in the correct position.
 - c. Inspect the rotor pins, hold the rotor up or turn it on its side and look into the rotor drive hole. you can see two metal pins in the hole (Figure 1 and 2).
2. Carefully lower the rotor straight down onto the drive spindle hub (Figure 3).
3. Make sure that you position the rotor is properly on the hub. (as shown in Figure 4).
4. Press the lid knob down on the rotor to fix it to the drive spindle hub. Turn the lid to the right (clockwise) until secure.

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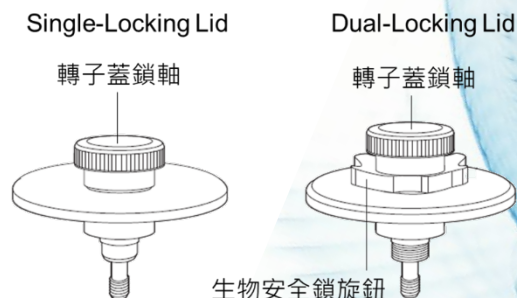
*Beckman Coulter 叫修專線: 0800-211-283

Confidential - Company Proprietary